


Traceable®

Thermohygrometer with Clock

Quick-Start Guide

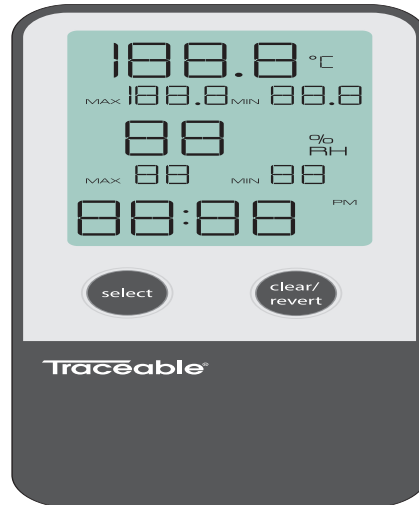

WARRANTY, SERVICE, OR RECALIBRATION

Traceable®
an antylia scientific company


Traceable

® Products are ISO 9001:2018 Quality-

Certified by DNV and ISO/IEC 17025:2017 accredited as a Calibration Laboratory
by A2LA.

Item No. 56000-10 Model No. 5660 ©2022 1065T6_M_92-5660-00 Rev. 0
09022022

DISPLAY DETAILS

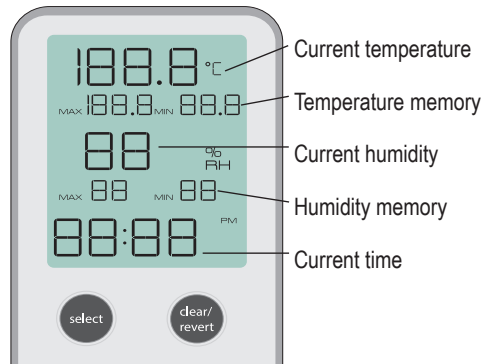


CLEARING INDIVIDUAL READINGS

1. Press select button until desired reading to be cleared is flashing.
2. Press and hold clear/revert button for 3 seconds to clear reading.

CLEARING ALL MEMORY READINGS

1. Press and hold clear/revert button for 3 seconds to clear all memory readings.

Traceable® Thermohygrometer
with Clock

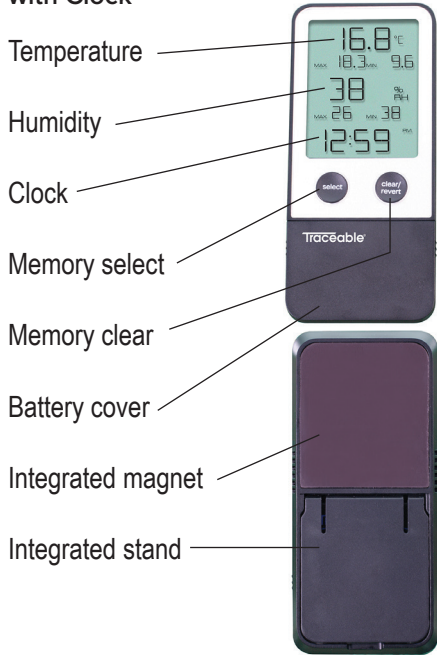


INITIAL SETUP


- 1 Remove battery cover and insert batteries.



- 2 Once batteries are installed, press O button to select setting.

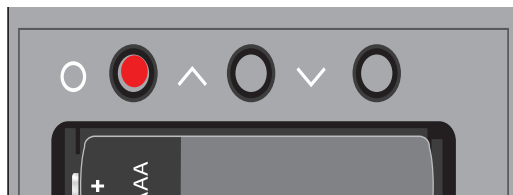

3 The first setting, Hours, will flash on display.

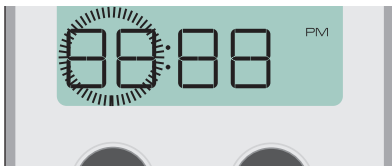

4 Press “^” and “v” buttons to adjust setting.



5 Once finished, press O button to save setting and proceed to Minutes, 12/24 Hour, etc.



6 Complete all settings to configure device:

- Hours
- Minutes

• 12 or 24 hour format clock

• Temperature units

7 Replace battery cover.

